

Hall D Detector Hall D Start Counter Checklist

Document ID: 466

Group: Hall D Detector

Subsystem: Hall D Start Counter

Author: Ito, Mark (marki)

Submitted By: Ito, Mark (marki)

Revision Number: 1

Revision Date: 2014-06-17 18:20

Revision Comment: first draft

1. check that bias voltages are on
2. check that low-voltage power is on
3. check that the start counter EPICS alarm handler is on
4. check that all EPICS alarms are clear
5. check that start counter temperatures are nominal
6. check that all counters are registering hits
 1. take data with the start counter cosmic trigger
 2. look the start counter monitoring histograms and check that they are nominal